**Méthode ESAO<sup>®</sup>**

***ESAO<sup>®</sup> Method***

**Ref :  
451 038**

**Français – p 1**

**English – p 17**

**Version : 7001**

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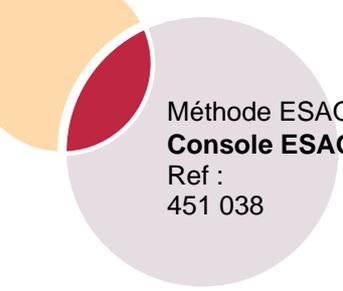
## 1 Introduction

The ESAO Primo is a computer-assisted experimentation console for laboratory use. To operate, it must be connected to the PC via a USB port. It has 4 input ports for VTT type sensors, and 4 direct differential  $\pm 20$  V input ports for use with analog output measuring apparatus, among other things.

It works with the *Atelier Scientifique* (Science Workshop) software (Compliss, Générés 5, Sérénis, Titration, Chronoméca, etc.).

To ensure a perfect supply of all sensors, the ESAO PRIMO console must be supplied using the mains adapter delivered with the console.

## 2 USB technology

Nowadays, conventional peripherals (mouse, printers, cameras, other photographic apparatus, etc.) are all equipped with a USB connection. This technology allows one to take advantage of a power and a use flexibility unlike any other before.

The USB port is found on all new computers. It is the next essential natural step in the development of your computer hardware.

The ESAO Primo console is compatible with most of your equipment and allows you to benefit from all of the advantages of USB technology (high speed, easy connection, etc.) by assuring the durability of your hardware with your future computer equipment.

## 3 Description

### Front side

## Back side

### 3.1 Sensor ports

There are four ports on the front side of the case for four separate VTT type sensors. All VTT type sensors will fit in these ports .

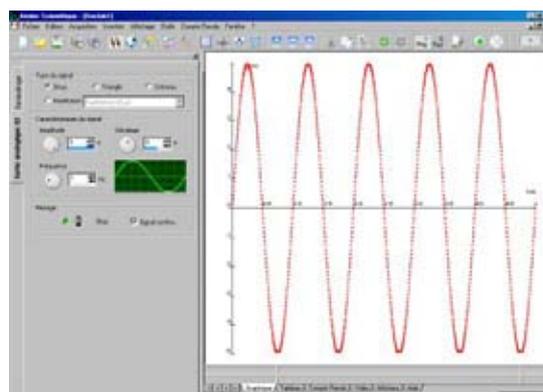
### 3.2 Direct analog input ports

The console is equipped with four  $\pm 20$  V differential analog input ports, which are accessible via 4 mm diam. safety sockets located on the back side of the console: a red socket for the + pole and a black one for the - pole. With these ports, the voltage of a current generated by an analog output measuring device, for example, can be measured directly. It is thus possible to measure a function generator signal. The signal from the analog output can be measured directly with the Générés 5 software. These input ports are protected against surges up to 250 V.

### 3.3 Analog output port

The ESAO Primo has a  $\pm 5$  V analog output port on the left side of the console that is accessible via 4 mm diam. banana safety sockets (yellow and black sockets).

This port can be accessed directly in the acquisition workshop of Générés 5. The signal generated can be measured directly on one of the ESAO console's direct differential input ports.

This output port can be used for certain applications that require the generation of a signal, such as the electrical activity of a crab nerve with the nerve-muscle software in SVT or an MPI application in Physics-Chemistry.

### 3.4 Synchronization input port

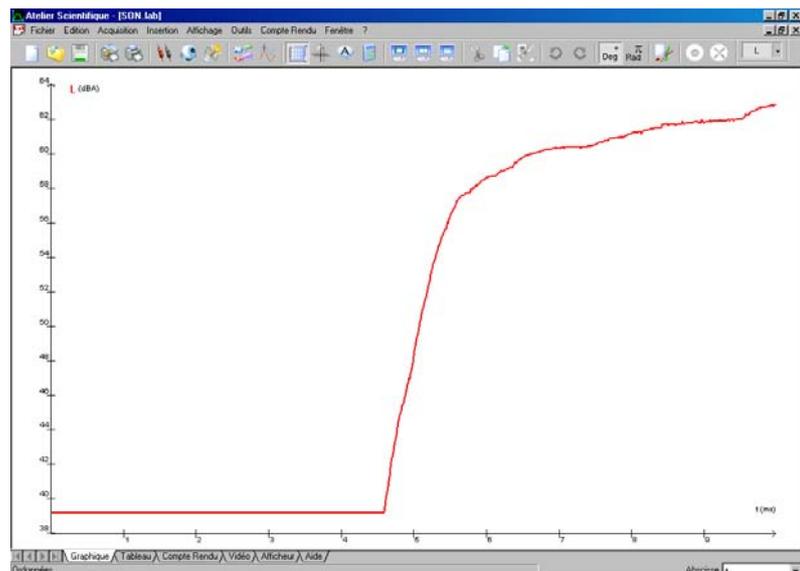
With this input port, you can launch an acquisition of a mechanically triggered event, such as one triggered by the Martoreflex in the myotatic reflex study in SVT or the study of sound in Physics-Chemistry.

This port is accessible via the red 4 mm diam. safety socket on the side of the console. The 0 V reference is provided via the adjacent 4 mm diam. black safety socket.

#### Example:

- Connect a sound sensor into one of the ports on the front side of the ESAO Primo console. This sensor may be used in either the signal mode or in the sound intensity mode in dB, as in the example below.
- Connect 2 safety cables with banana plugs to the synchronization input port sockets (1 in the red socket and 1 in the black socket).
- Attach the metal rods to the ends of the banana cables in order to create a clap function.
- Set the software parameters for a positive edge-triggered acquisition of 10 ms and 1000 points.
- Launch the acquisition by clicking on the green light, which is now waiting for the synchronization.
- Striking the 2 rods together creates a short circuit that generates a reduction in voltage from 0 to 5 V at the synchronization input port, indicating the start of the measurement.

#### Example of the results:

## 4 Driver installation

It is necessary to install a driver to operate the console, which is found on the CD ROM shipped with the console. The driver only has to be installed with the initial use.

### 4.1 In Windows 2000

1. Make sure that the cable is not plugged into the USB port.
2. Open a session in the **administrator** mode (i.e., use an account with administrative rights).
3. Plug the cable into a USB port on the computer.
4. Wait a few seconds until the PC detects the new peripheral. A window will open automatically once it is detected.

5. Insert the CD ROM in the disk drive if you have not already done so.
6. Click the *Next* button to continue the installation.

7. Select the search for a suitable driver for my device (recommended) option and click Next.

The system will search for the right driver.

8. The peripheral and the driver location are now identified. Click Next.
9. Windows will copy the files (this may take several minutes) and will indicate that it has completed the installation.

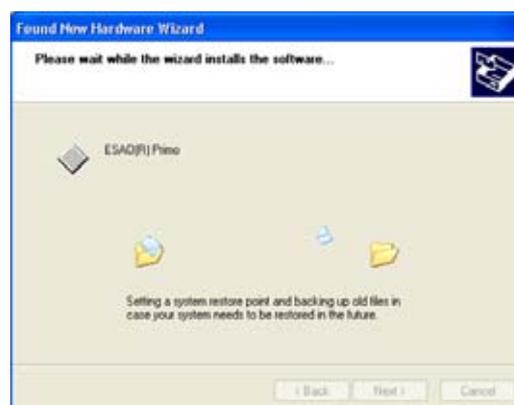
Click Finish.

## 4.2 In Windows XP

1. Make sure that the cable is not plugged into the USB port.  
Open a session in the **administrator** mode (i.e., an account with administrative rights).
3. Insert the CD ROM in the disk drive if you have not already done so. Click on the No, not at this time option and then click the Next button.
4. If nothing happens in 30 seconds after inserting the CD ROM, select Install the software automatically (recommended) and click Next.

5. Windows will search for drivers for the USB cable. If an intermediate ("Hardware installation") window opens, you must click the *Continue* button.
6. Windows will copy the files (this may take several minutes).

7. Windows will indicate that it is finished when the installation is complete. Click Finish.

## 5 Software compatibility

The ESAO Primo is compatible with the software in the Science Workshop collection:

Généris 5, Sérénis, Compliss, Actocard, Actospir, Actoresp, Actotensio, Nerf-Muscle, Actohom, Chronoméca, Titrage.

### **It will not work with the 16 bit SVT software**

Respi, QR, Methom, Respihom, Neuro, Cardio, Spirom, Reacell, Cell-C, Phyeff, Tensio, Enzymo, Refmyo, Winrad

### **It will not work with the "junior software"**

Echanges junior, Souffle junior, Animo-junior, Cardio-junior for the SVT and Watt-Ohm junior, pH junior, Calor junior

## 6 Operating

The VTT type sensors are recognized automatically by the console and the Science Workshop software.

## 7 Physics-Chemistry acquisition example

### 7.1 Data acquisition with the Compliss software: Ohm's law

Start Compliss or Généris 5. Drag the voltmeter  to one of the lines of the Y axis and designate the voltage U. The sensor selector switch should be on continuous mode.

**You can use a direct differential input port on the ESAO Primo console. If you do so, there is no need to use the VTT voltmeter sensor. In this**

**case, drag the  icon symbolizing the direct input port to one of the lines on the Y axis.**

Drag the amperemeter  to the X axis. Designate the current I. Set the graph interval between 0 and 150 mA.

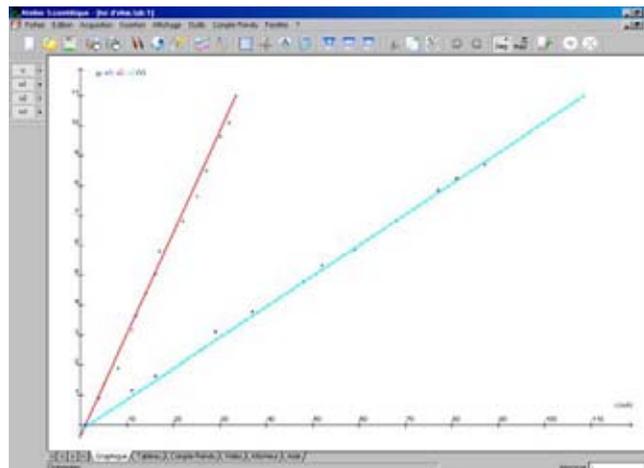
Click on the Validation tab. Choose automatic with 1% spread.

Click on  to launch the acquisition and then vary the voltage.

Click on  to stop the acquisition.  
 File, save as... Select the file folder and then name it **Ohm.lab**.

## 7.2 Results

## 8 SVT acquisition experiments

### 8.1 Data acquisition with the Compliss software: Gas exchanges in plants

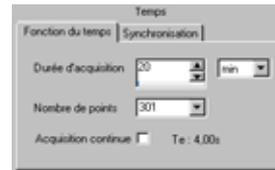
Start Compliss or Sérénis.

Drag the oxymeter sensor  to one of the lines on the Y axis and then do the same with the luxmeter.

If necessary, calibrate the probe.


Drag the time icon  Temps to the X axis.

Set the parameters to 20 minutes and 301 points :

Place the experimental device in a dark place and launch the acquisition by clicking on . Wait a few minutes, and then expose the device to full light with a cold lamp. Repeat the operation several times.

## 8.2 Results
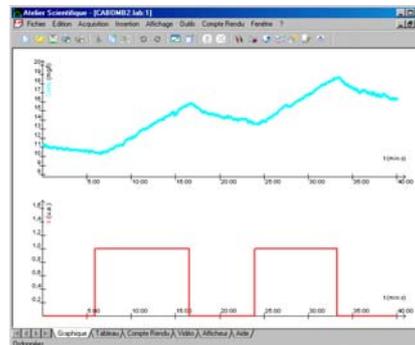
## 9 Technical specifications

### 9.1 Analog input ports

- Sensor input port:
  - Resolution: 12 bits
  - Accuracy: +/- 0.2%
  - Conversion time: 10 μs
  - Maximum effective conversion rate: **100/n kHz on n ways**
- Direct input ports:
  - Scale: +/-20V
  - Resolution: 12 bits
  - Accuracy: +/- 0.2 %
  - Conversion time: 10 μs
  - Maximum effective conversion rate: **100/n kHz on n ways**

### 9.2 Analog output port

- 1 analog output port** accessible via the 4 mm yellow banana sockets:
- Scale: +/-5 V
  - Resolution: 12 bits
  - Minimum conversion time: 10 μs
  - Maximum calibration frequency: 100 kHz

### **9.3 External SYNCHRO input port**

**1 synchronization input port** accessible via the red 4 mm banana socket:

- Level: 0 – 5V
- Triggering: Positive or negative edge

### **9.4 Power supply**

The power supply is ensured by the mains block delivered with the console :  
100 à 240 V/~ /9 V (regulated  $\pm 5\%$ ;  $\geq 1.3$  A.

### **9.5 Operating conditions**

#### **9.5.1 Temperature**

- Operating : from 0 to 40°C max.
- Avoid prolonged exposure to sunlight.

#### **9.5.2 Electromagnetic compatibility**

The measurements may be momentarily distorted under certain unusual electromagnetic environmental conditions. This in no way affects the accuracy of the apparatus.

## **10 After-sales service**

The apparatus is covered by a 2 year warranty. In the event of defects, it should be returned to our shops.

For all repairs, adjustments, or replacement parts, please contact:

**JEULIN TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**+33 (0)2 32 29 40 50**

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

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We declare under our own responsibility, that the following products :

**ESAO® Primo ref. : 451 038 associated to the AC adaptor and to his sensing adaptors**

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