# METHODE VTT

Capteur VTT

## *VTT Method*

*VTT sensor*

Ref :  
472052

Français – p 1

English – p 5

Version : 6007

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***VTT ammeter sensor 2.5A / 10A***

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**NOTES**

## 1 General information

This sensor allows to measure intensities of current with a VTT system or ESAO Primo. Measurements are carried out by a differential entry with safety banana plugs.

## 2 Use

### 2.1 Interfaces

This sensor operates with the graph console or the alphanumeric VTT console or ESAO Primo.

### 2.2 Connecting the sensor to the interface

The sensor is connected by means of a connector located at the back of the box:

- Place the sensor in line with the selected measurement channel.
- Engage the sensor on the console guide,
- follow the guidance system of the VTT console ,
- Engage it fully.

### 2.3 Type of measurement

The sensor measures the instantaneous value of the current passing through.

### 2.4 Scale

The user has 2 scales of measurement:                   -2.5 /+ 2.5 A  
  -10 A /+ 10 A

The scale is selected by means of the switch on the box.

### 2.5 Startup with VTT

- connect the sensor,
- light up the VTT console.

If the sensor is not detected, update your VTT console connected to the PC with the CD via the ESPACE-VTT software.

- connect the console to your computer

- open Espace VTT by clicking on the icon  displayed on your workstation

- click on install:

- install the Biosflash software by clicking on open.

- The VTT console is then automatically updated
- disconnect your VTTconsole : the sensor is then detected.

## 2.6 Startup with ESAO Primo

The sensor must be recognized by the interface.

- connect the console to your computer,
- connect the sensor on the console,
- start your acquisition software.

## 2.7 Technical specifications

- Measuring range: from - 2,5 A/+ 2,5 A  
from -10 A /+ 10 A
- Resolution: 5 mA  
10 mA
- Precision: 1% to+/- 15 mA 2.5 A caliber,  
1% to+/- 60 mA 10 A caliber
- Band-width: 0-10 Khz
- Resistance of entry: 0.01 Ohm
- Protection: Fuse
- Tension common mode: 35 V

### **3 After-Sales Service**

This material is under a two year warranty and should be returned to our stores in the event of any defects.

For any repairs, adjustments or spare parts, please contact:

**JEULIN - TECHNICAL SUPPORT**  
**Rue Jacques Monod**  
**BP 1900**  
**27 019 EVREUX CEDEX FRANCE**  
**+33 (0)2 32 29 40 50**

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediately to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

**Free service \* :**  
**+ 33 (0)2 32 29 40 50**

*\* Call cost not included*