**Méthode VTT<sup>®</sup>**

***VTT<sup>®</sup> method***

**Ref :  
472 021**

**Français – p 1**

**English – p 5**

**Version : 6005**

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***VTT<sup>®</sup> sound level sensor***

## 1. Description

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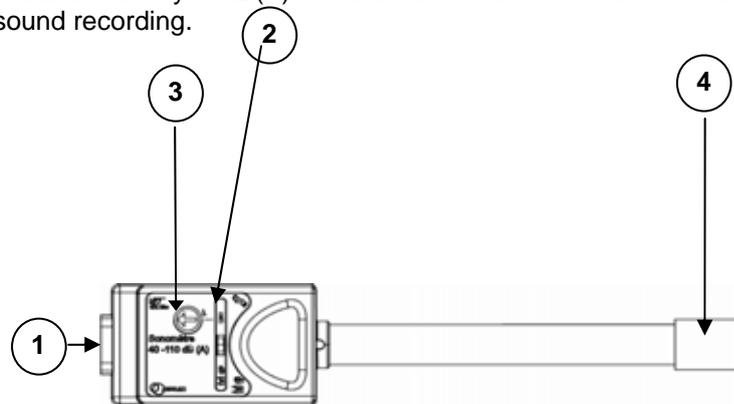
**NOTES**

## 1. Description

The VTT sound level sensor enables the measurement of an audible signal or a sound intensity in dB(A). It includes an electret-condenser microphone for sound recording.

1. Connection to the console
2. Function selector switch: sound or sound-level metre.
3. Potentiometer for adjusting the audio amplitude for the sound function.
4. Protective cover for the electret-condenser microphone.

## 2. Use

### 2.1 Interfaces

This sensor works with the VTT<sup>®</sup> console or ESAO Primo.

### 2.2 Connecting the sensor to the VTT<sup>®</sup> console

The sensor is connected using the connector located at the back of the box of the VTT console or the front face of the ESAO Primo:

- Place the sensor in line with the selected measurement channel.
- Engage the sensor on the console guide.
- Follow the guidance system of the VTT<sup>®</sup> console.
- Engage it fully.

It can then be used with VTT:

- In stand-alone mode

or

Connected to a computer (an acquisition software such as Generis is then required).

## 2.3 Commissioning with the VTT console

- Connect the sensor.
- Switch on the VTT® console.

If the sensor is not detected, update your VTT console connected to the PC using the floppy disc via the Espace VTT software.

- Connect the console to the PC,
- Open the Espace VTT by clicking on the icon  present in “My computer”,
- Click on Install ,
- Install the Biosflash software.

## 2.4 Commissioning with the ESAO Primo

The sensor is automatically detected by the sensor.

- Connect the console to your PC,
- Connect the sensor to the console,
- Start your acquisition software.

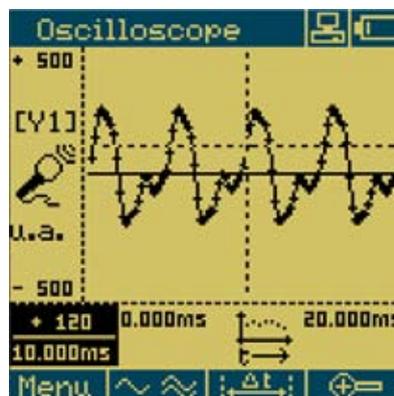
## 3. Technical characteristics

<b>Functions:</b>	By switch (2): <ul style="list-style-type: none"><li>• Sound, +/- 500 a.u. (arbitrary unit)</li></ul> Resolution 1 a.u. <ul style="list-style-type: none"><li>• Sound level meter, 40 to 110 dB(A)</li><li>• Accuracy +/- (5% + 0.21dB(A))</li><li>• Resolution 0.07 dB(A)</li></ul>
<b>Adjustment</b>	Adjustment of the amplitude of the audible signal (3). The adjustment enables the adaptation of the amplification to the sound intensity received.

**Note:** do not immerse the microphone in water.

## 4. Experimentation

It is possible to display and record the shape of any audible signal in sound mode:

In sound level meter mode, it is also possible to measure and record sound intensities.


## 5. After-sales service

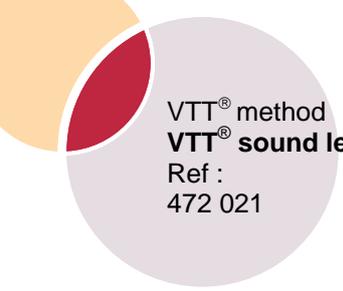The device is under a 2-year guarantee, it must be sent back to our workshops.

For any repairs, adjustments or spare parts, please contact:

**JEULIN – TECHNICAL SUPPORT**  
Rue Jacques Monod  
BP 1900  
27 019 EVREUX CEDEX FRANCE  
+33 (0)2 32 29 40 50

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SA capital 3 233 762 € - Siren R.C.S. B 387 901 044 - Siret 387 901 04400017

## Direct connection for technical support

A team of experts at your disposal from Monday to Friday (opening hours)

- You're looking for technical information ?
- You wish advice for use ?
- You need an urgent diagnosis ?

We take in charge your request immediatly to provide you with the right answers regarding your activity field : Biology, Physics, Chemistry, Technology .

**Free service \* :**  
**+ 33 (0)2 32 29 40 50**

*\* Call cost not included*