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REF. 472 004

I - GENERALITES

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REF. 472 004

I - OVERVIEW

This sensor enables measurement of temperature from -20 to 120 °C with the VTT system.

The probe used is a 316L stainless steel probe (pharmaceutical quality) of PT1000 type.

II - USE

1. VTT system

The sensor works with the **VTT console**: Ref. 477 000

2. Connecting the sensor to the VTT console

The sensor connects by means of the connector located at the back of the casing:

- introduce the sensor aligned with the measurement channel chosen.
- locate the sensor in the **console**'s guide slots.
- follow the VTT **console**'s guide slot system.
- locate fully home.

3. Scale

The user has a measurement scale from -20 to 120 °C.

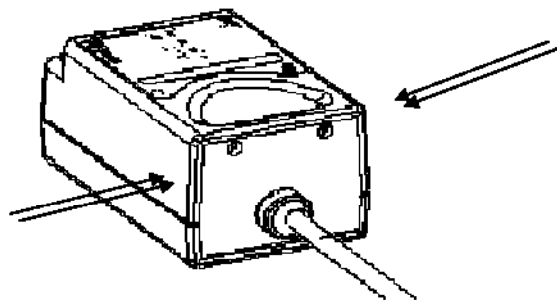
4. Use

- connect the sensor.
- switch on the VTT console.

5. Calibrating the probe

In the event of deviation in measurements, proceed as follows:

- prepare a recipient, R0, containing a bath of melting ice (at 0°C).
- prepare another recipient, R100, containing pure, boiling water (at 100 °C).
- locate the potentiometers' zeroing access orifices on the front of the sensor (P100 at left and P0 at right).
- connect the sensor.
- switch on the VTT console.
- use the screwdriver supplied with the console to make the following adjustments:
 - 1 - Plunge the probe into R0.
 - Wait for the measurement to stabilise.
 - Set P0 to display 0 °C.
 - 2 - Plunge the probe into R100.
 - Wait for the measurement to stabilise.
 - Set P100 to display 100 °C.

Repeat points 1 and 2 several times to simultaneously get measurements that are accurate in line with the values stated in paragraph III.

III - TECHNICAL FEATURES

- | | |
|---------------------------------|---------------------|
| - Length: | 135 mm |
| - Diameter: | 6 mm |
| - Measurement range: | from -20 to 120 °C. |
| - Resolution: | 0.14 °C. |
| - Accuracy: | ± 2 °C. |
| - Response time in water: | about 15 s. |
| - Maximum exposure temperature: | -50°C./+200°C |

IV - CARE AND MAINTENANCE

For any repairs, adjustments or spare parts, contact:

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